

AgriSpec® User Manual



ASD Document 600542 Rev. E

© 2008 by ASD, Inc.
www.asdi.com



Trademark Information

ASD, Inc.

2555 55th Street Suite 100
Boulder, CO 80301 USA
Phone: (303) 444-6522
www.asdi.com

FieldSpec®, LabSpec®, QualitySpec®, RxSpec®, ViewSpec™, Indico™, AgriSpec®, RS³™, and TerraSpec® are registered and unregistered trademarks, and the intellectual property of ASD, Inc. (ASD), Boulder, Colorado, USA. All trademarks used or displayed in this document are trademarks of ASD, its affiliates, or third party owners. Unauthorized use of these trademarks is illegal and punishable by law. Nothing contained in this document is to be construed as granting, by implication, estoppel, or otherwise, any license or right of use of any such trademark without the prior and express written permission of ASD, or such third party owner.

This document contains proprietary information protected by copyright law and may not be reproduced in any manner without the express written approval of ASD, Inc.

The information and specifications contained in this manual are subject to change without notice. ASD, Inc. shall not be held liable for technical, editorial omissions, or errors made herein; nor for incidental or consequential damages resulting from furnishing, performance, or use of this material.

Technical Support

If you have any questions or concerns, please contact ASD, Inc. by phone, fax, or email:

Phone: 303-444-6522 X-144
Fax: 303-444-6825
email: support@asdi.com
Web: www.asdi.com
ftp: [ftp.asdi.com](ftp://ftp.asdi.com)

Technical support is committed to providing you with a timely response to your questions. We will work with you to provide solutions to your applications. Technical support is available to answer your questions Monday thru Friday, 8 am to 5 pm Mountain Standard Time. We will happily respond to your e-mail queries as well.

Table of Contents

	Trademark Information	ii
	Technical Support	ii
	Declaration of Conformity	vii
Chapter 1	Introduction	1
	Hardware Specifications	3
	Physical Specifications	3
	Power Input and Output	3
	AgriSpec Wavelength Configuration	4
	Interface Requirement	5
	Standard Accessories	5
	Computer Requirements	6
	Software Requirement	6
	Theory of Operation	7
Chapter 2	Setup	11
	Unpacking the Instrument	12
	Power Supply	13
	Ethernet Cables	15
	Accessory Power Port	16
	Remote Trigger	17
Chapter 3	Usage and Maintenance	19
	Light Sources	19
	Access and Ventilation	19
	Cleaning	20
	Annual Maintenance	20
	Returning Instrument to ASD for Service	20
Chapter 4	Fiber Optic Interface	21
	Fiber Optics Cables	21
	Fiber Optic Cable Interface	22
	Matching Fiber Bundles	24
	LED Check for Fiber Optic Cable	26

Chapter 5	Batteries	31
	Battery Care	31
	Battery Charger	32
	Application Battery Indicator	33
Chapter 6	Field Measurements	35
	Illumination	36
	Vegetation	36
	Rocks, Soils, and Man-Made Materials	37
	Suggested Set-up	37
	RS3 Application	37
	Indico Application	38
	White Reference	39
	Spectralon Reflectance Data	39
	Maintaining Spectralon References	41
	White Reference Procedures	42
Appendix A	Troubleshooting	43
	Common Communication Fixes	43
	Does Not Connect to the Spectrometer	43
	Does Not Connect Wireless	45
	Does Not Connect to the Access Point	46
	Spectrometer Loses its Wireless Connection	47
	Wireless Interface is Disabled in the Netcfg Application	47
Appendix B	Reference Information	49
	Frequently Asked Questions (FAQ)	49
	What is a spectrometer?	49
	What are the differences between the ASD spectrometers?	50
	What is remote sensing?	50
	How often do I need to optimize?	51
	How often do I need a (baseline) White Reference?	51
	What spectrum average (or sample count) should I use?	51
	What if I want to keep the same settings all day?	52
	What are the units of radiance?	52
	Can I post-process my data?	52
	Why do I see oscillations (sine wave) in my data?	52
	What are these upward or downward spikes in VNIR data?	52
	What are these steps in my data?	53

What can cause more noise in my data from last time?	53
Why does the VNIR drop to zero after a Dark Current collection?	54
How do I collect a reference with a spot size larger than the Spectralon panel?	54
How do I set up GPS?	54
What does a broken fiber mean?	55
How long is the battery life?	56
What type of Ethernet cable can I use for the static IP configuration? . .	56
Can I install additional software on the instrument controller?	57
Why does the software seem to do unexpected things?	57
How long does it take for the AgriSpec spectrometer to warm up?	57
Where is my serial number?	57
Data Conversion	58
Gain and Offset	58
Optimization	59
Spectrum Averaging (or Sample Count)	59
Temperature Effects	59
Technical Support	60
W.E.E.E. Compliance	60
Index	61

Declaration of Conformity

According to IEC guide 22 and EN45014

Manufacturers Name: ASD Incorporated.

Manufacturers Address: 2555 55th Street, Suite 100, Boulder, Colorado 80301.

Declares that the product

Product Name: AgriSpec™

Product Numbers	Description	Range (nm)
A100900	Full Range (VNIR, SWIR1, & SWIR2)	350-2500
A100910	SWIR2 only	1800-2500
A100920	SWIR1 and SWIR2	1000-2500
A100930	SWIR1 only	1000-1800
A100940	VNIR only	350-1050
A100950	VNIR and SWIR1	350-1800

Conforms to the following EU Directives:

Safety: Low Voltage Directive, 72/23/EEC, as amended by 93/68/EEC

EMC: Electromagnetic Compatibility Directive, 89/336/EEC, as amended by 93/68/EEC

Supplementary Information:

The product complies with the requirements of the following Harmonized Product Standards and carries the CE-Marking accordingly:


EN 61010-1: 2001

Safety Requirements for Electrical Equipment for Measurement, Control and Laboratory use

EN 61326: 2002

Class A, Electrical Equipment for Measurement, control and Laboratory use – EMC requirements

Signature:

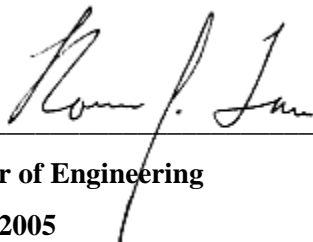

Title: Director of Engineering

Date: 11-02-2005

For compliance Information ONLY, contact:

European Contact: Your local ASD Representative.

USA Contact: Product Regulations Manager, ASD Inc.

2555 55th Street, Suite 100, Boulder, Colorado 80301.

Phone: (303)-444-6522.

Symbols - Definitions



CAUTION: *Risk of danger.* This is a personal danger warning. Documentation must be consulted in all cases where this symbol is marked. Failure to acknowledge these warnings could result in personal injury to the user.



CAUTION: *Risk of Electric Shock.* This is a personal danger warning. Documentation must be consulted in all cases where this symbol is marked. Failure to acknowledge these warnings could result in personal injury to the user.



CAUTION: *Hot Surface.* This is a personal danger warning. Documentation must be consulted in all cases where this symbol is marked. Failure to acknowledge these warnings could result in personal injury to the user.



Recycle. Items with this symbol indicate that the item should be recycled and not disposed of as general waste.

Warnings and cautions are placed throughout this manual for the convenience of the reader. However, the absence of warnings and cautions do not preclude the use of proper caution and handling. Usual precautions are recommended to be taken at all times, either written or otherwise, to avoid personal injury or damage to ASD equipment.

The Power Supply is intended for Indoor Use Only

When using an electrical appliance, basic safety precautions should always be followed.

Danger- To reduce the risk of electric shock read all instructions before using the power supply. A power supply should never be left unattended when plugged in. Always unplug the power supply from the mains socket immediately after using.

Warning- To reduce risk of burns, fire, electric shock or injury to persons or animals:

1. Use this power supply only for its intended use as described below.
2. Do not use outdoors.
3. Do not allow to be used as a toy. Pay close attention when this power supply is used by, or near to, children.
4. Use only attachments recommended by the manufacturer.
5. Never operate the power supply if it has a damaged cord or plug, if it has been dropped or damaged or if it has fallen into water. In such cases return the power supply to an authorized dealer or service center for examination or repair.
6. Never drop or insert an object into any openings.
7. Do not operate where aerosol (spray) products are being used or where oxygen is being administered.
8. The power supply should be used near to a convenient and easily accessible mains socket.

Notes:

Chapter 1 Introduction

The AgriSpec spectrometer is a general-purpose instrument that measures the visible near-infrared (VNIR) and short-wave infrared (SWIR) spectra for agriculture remote sensing and analysis.

The AgriSpec spectrometer is useful in many application areas, but in particular agriculture and the field analysis of vegetation and soils. The resolution is ideally suited for performing reflectance and transmittance spectral measurements of a wide range of samples.

The ASD Accessories Guide contains an array of sampling accessories for the AgriSpec spectrometer, such as the plant probe and leaf clip, that extend the application areas.

The AgriSpec spectrometer is compact and field portable. It has a spectral range of 350-2500 nm and rapid data collection (10 spectra per second).

While the most highly regarded features are performance and field-portability, these instruments perform competitively in the laboratory as well.

Features and Advantages

Accurate and Precise - High signal-to-noise ratio and superior repeatability of results for better discrimination and analysis of materials.

Rugged and Transportable - Resistant to vibration and changes in temperature or humidity.

Fast - 10 spectra per second data collection for the entire 350-2500 nm range.

- » 0.1 second scanning time provides an accurate average every second
- » Users can quickly scan several areas when analyzing bulk samples.
- » Rapid analysis allows measurement of large numbers of samples.

Compatible - Works with off the shelf chemometrics software (Grams™ PLS plus/IQ and Unscrambler®).

1.1 Hardware Specifications

This section contains details about the AgriSpec spectrometer.

1.1.1 Physical Specifications

Height	12.7 cm	5 inches
Width	35.6 cm	14.5 inches
Depth	29.2 cm	11.5 inches
Weight	5.6 kg	12 lbs

- Durable satin powder coat finish with green urethane end-caps and handles
- All vital components are in a dust-proof enclosure and EMI sealed
- Adjustable fiber optic to instrument spacer

1.1.2 Power Input and Output

AC Power Supply Type	Auto ranging, Switching, SELV
AC Input	90-240 VAC, 50/60 Hz
DC Input	+12 VDC, 50 Watt
Accessory Power Port (front of instrument)	Output, +12 VDC, 19.2 W (max).

The near infrared (NIR) is also called the short-wave infrared (SWIR).

Instrument Configuration	Power Rating (Maximum)
350-1050 nm	12 VDC, 45 Watt
350-1800 nm	12 VDC, 50 Watt
1000-1800 nm	12 VDC, 50 Watt
1800-2500 nm	12 VDC, 50 Watt
1000-2500 nm	12 VDC, 60 Watt
350-2500 nm	12 VDC, 60 Watt

1.1.3 AgriSpec Wavelength Configuration

Wavelength Name	Wavelength Range
VNIR-SWIR1-SWIR2	350-2500 nm
VNIR only	350-1050 nm
VNIR-SWIR1	350-1800 nm
SWIR1 only	1000-1800 nm
SWIR1-SWIR2	1000-2500 nm
SWIR2 only	1800-2500 nm

The spectral resolution is:

- 3 nm (Full-Width-Half-Maximum) at 700 nm.
- 10 nm (Full-Width-Half-Maximum) at 1400 nm.
- 10 nm (Full-Width-Half-Maximum) at 2100 nm.

The sampling interval is:

- 1.4 nm for the spectral region 350-1000 nm.
- 2 nm for the spectral region 1000-2500 nm.

The AgriSpec spectrometer may be configured to have three separate holographic diffraction gratings with three separate detectors. Each detector is also covered with the appropriate order separation filters to eliminate second and higher order light.

- VNIR: 512 element silicon photo-diode array for the spectral region 350-1000 nm.

- SWIR1: graded index, TE-cooled, extended range, InGaAs, photo-diode for the spectral region 1000 nm to 1800 nm.
- SWIR2: graded index, TE-cooled, extended range, InGaAs, photo-diode for the spectral region 1800 nm to 2500 nm.

1.1.4 Interface Requirement

The AgriSpec spectrometer has 10/100 Base T Ethernet port. The spectrometer with an Ethernet cross-over cable can be directly connected to the Ethernet port provided on many laptop computers.

The AgriSpec spectrometer can communicate with the instrument controller using an optional 802.11b wireless card. If the instrument controller is supplied by the customer, it must be 802.11b compatible.

The AgriSpec spectrometer uses industry standard components. It has the same general connection speed and distance capabilities as other 802.11 lb wireless devices.

For the best performance, ensure that there is a clear line-of-sight between the spectrometer and instrument controller. Obstructions or radio frequency interference in close proximity can cause a loss of communication or significantly shorten the communication range.

Note: For information about configuration of the Ethernet card and cable, please refer to the *Installation Manual*.

1.1.5 Standard Accessories

- AgriSpec™ Spectrometer
- AgriSpec® User Manual (PDF on Flash Drive)
- Instrument Controller: ASD-approved laptop computer
- RS³ Software Pack
- Indico Software Pack
- 1.5 meter FR Low OH Fiber Optic Jumper Cable
- Power Supply 12 V
- Battery Belt, Gel Cell (12 V, 9 A hour) with Charger and 6 ft. Battery-to-Instrument Power Cable
- Power Adapter with 6 m (20 ft) Cable for Automobile Cigarette Lighter
When using this adapter, turn off the engine to the vehicle, electromagnetic fields from the engine can generate noise in the acquired signals.
- 1.5 meter Instrument-to-Probe Accessory Power Cable

- Remote Trigger with LEDs and Velcro Pads
- RJ45 CAT 5e UTP Ethernet, Shielded, Crossover Cable
- 64 MB (min.) Flash Drive Memory USB Flash Drive
- Mylar Wavelength Reference
- Magnifier for Fiber Checker
- Shipping Trunk: Durable, air tight, wheeled, fold down handle, and foam filled with cutouts to protect the instrument and many accessories

1.2 Computer Requirements

The *instrument controller* is a computer that manages the AgriSpec spectrometer, stores data, and processes the results.

The standard AgriSpec spectrometer includes a laptop computer, or the *instrument controller*.

The minimum requirements for the instrument controller are:

- 1.2 GHz Pentium or better notebook or PC-w/monitor
- 256 MB RAM or more
- 20 GB of free disk space
- 1024 x 768 or better graphics resolution
- 24-bit color or better - 32-bit recommended
- Ethernet port: 10/100 Base T Ethernet interface
- (Optional) Ethernet wireless (WiFi) adaptor: PCMCIA, USB, or built-in that is compatible with the 802.11b standard
- (Optional) Serial communications port (or USB port) for GPS receiver (Only if a GPS receiver will be used)
- (Optional) NMEA compatible GPS receiver

1.3 Software Requirement

The instrument controller requires the following software:

- Either RS³ Software from ASD
- Or Indico Software from ASD
- Microsoft Windows® 95/98/NT/2000/ME/XP Operating System

- Microsoft Internet Explorer 6.0 or better.

Users need a basic understanding of the Microsoft Windows operating system including software installation.

International customers using non-English versions of Windows must alter the **Regional Settings** under **Start->Settings->Control Panel**. The default language must be set to English (United States) in order for the software to be registered and operate correctly. The numbering format must also be set to English.

1.4 Theory of Operation

The AgriSpec spectrometer measures the optical energy that is reflected by, absorbed into, or transmitted through a sample. *Optical energy* refers to a wavelength range that is greater than just the visible wavelengths, and is sometimes called *electromagnetic radiation* or *optical radiation*.

With accessories, various set-ups, and built-in processing of the optical energy signal, the AgriSpec spectrometer can measure:

- spectral reflectance,
- spectral transmittance, and
- spectral absorbance.

Fiber Optic Collection of Reflected/Transmitted Light

Optical energy is collected through a bundle of specially formulated optical fibers, precisely cut, polished, and sealed for extremely efficient energy collection. The fibers themselves are of low OH composition providing the maximum transmission available across the wavelength range of the instrument.

Inside the AgriSpec™ Spectrometer

The fiber cable delivers the collected optical energy into the instrument, where it is projected onto a holographic diffraction grating. The grating separates and reflects the wavelength components for independent measurement by the detectors.

Visible/Near-Infrared (VNIR)

The Visible/Near-Infrared (VNIR: 350-1000 nm wavelength) portion of the spectrum is measured by a 512-channel silicon photodiode array overlaid with an order separation filter. Each channel (or detector) is geometrically positioned to receive light within a narrow (1.4 nm) range. The VNIR spectrometer has a spectral resolution (full-width half maximum of a single emission line) of approximately 3 nm at around 700 nm.

Each detector converts incident photons into electrons. This photocurrent is continually converted to a voltage and is then periodically digitized by a 16-bit analog-to-digital (A/D) converter. This digitized spectral data is then transmitted to the instrument controller for further processing and analysis by the controlling software.

The 512-channel array permits the entire VNIR spectrum to be scanned in parallel at 1.4 nm wavelength intervals. A single sample can be acquired in as little as 17 ms.

Short-Wave Infrared (SWIR)

The Near-Infrared (NIR), also called Short-Wave Infrared (SWIR), portion of the spectrum is acquired with two scanning spectrometers:

- SWIR1 for the wavelength range of 1000 nm to 1800 nm
- SWIR2 for the wavelength range of 1800 nm to 2500 nm

The SWIR scanning spectrometers only have one detector for SWIR1 and another for SWIR2. This is different from the VNIR spectrometer, which has an array of 512 detectors. Thus, SWIR spectrometers collect wavelength information sequentially rather than in parallel.

Each SWIR spectrometer consists of a concave holographic grating and a single thermo-electrically cooled Indium Gallium Arsenide (InGaAs) detector. The gratings are mounted about a common shaft which oscillates back and forth through a 15 degree swing. As the grating moves, it exposes the SWIR1 and SWIR2 detectors to different wavelengths of optical energy. Each SWIR spectrometer has ~600 channels, or ~2 nm sampling interval per SWIR channel. The spectrometer firmware automatically compensates for the overlap in wavelength intervals.

Like the VNIR detectors, the SWIR1 and SWIR2 detectors convert incident photons into electrons. This photocurrent is continually converted to a voltage and is then periodically digitized by a 16-bit analog-to-digital (A/D) converter. This digitized spectral data is then transmitted to the instrument controller for further processing and analysis by the controlling software.

The grating is physically oscillating with a period of 200 ms. It performs a forward scan and a backward scan, resulting in 100 ms per scan. This is the minimum time required for any SWIR samples, or full-range samples.

Communicating with the Instrument Controller (Computer)

The AgriSpec spectrometer communicates with the instrument controller using a cross-over Ethernet cable or optional Wireless Ethernet (WiFi) interface. The amount of data that is sent depends on the configuration of the spectrometer.

- A single sample of VNIR is ~1024 bytes
- A single sample of SWIR1 is ~2400 bytes
- A single sample of SWIR2 is ~2400 bytes

A full-range AgriSpec spectrometer can create packet data sizes over 5 KB. Other single or dual range configurations of the spectrometer create smaller data packets (adding the packet sizes as above.)

When features for spectrum averaging (or sample count) are turned on within the application software on the instrument controller, the averaging is performed at the spectrometer.

Dark Current Measurement

Dark Current (DC) refers to current generated within a detector in the absence of any external photons. DC is the amount of electrical current that is inherent in the spectrometer's detectors and other electrical components and is additive to the signal generated by the measured external optical radiation.

Note: *Noise* is the uncertainty in a given measurement, one channel at a time. Noise by definition is random. Noise can be reduced by using more samples and averaging the signal. *Dark Current* is different from noise, because it is relatively stable and can be characterized.

This manual uses DC to refer to all systematic contributions to the detector signal. DC is a property of the detector and the associated electronics (not the light source). DC varies with temperature. In the VNIR region, DC also varies with integration time.

Whenever DC is taken, a mechanical shutter is used to block off the entrance slit of the VNIR spectrometer so the signal can be measured. This signal is subtracted from each subsequent spectrum until another DC is taken. The SWIR spectrometers take and subtract DC on every scan.

The DC measurement can be updated at any time, but should be updated more frequently in the beginning of a given session while the instrument warms up.

The VNIR spectrometer is fitted with a unique software and hardware combination feature called *Driftlock*. Driftlock corrects for DC changes over time. It automatically updates DC for every measurement by looking at a series of masked pixels at the front portion of the VNIR array. The Driftlock feature corrects for the majority of DC variation over time.

White Reference

A material with approximately 100% reflectance across the entire spectrum is called a white *reference panel* or white *reference standard*.

The raw measurement made by the spectrometer is influenced by both the sample and the light source. An independent measure of the light source illumination on a reference of known reflectance is required to calculate the reflectance of the sample. The use of a white reference standard with near 100% reflectance simplifies this calculation.

Reflectance and transmittance are inherent properties of all materials and are independent of the light source.

Reflectance is the ratio of energy reflected from a sample to the energy incident on the sample. *Spectral Reflectance* is the reflectance as a function of wavelength.

Transmittance is the ratio of the radiant energy transmitted through a sample to the radiant energy incident on the surface of the sample. *Spectral Transmittance* is the transmittance as a function of wavelength.

Relative reflectance is computed by dividing the energy reflected from the sample by the energy reflected off a white reference panel or standard.

Accessories for Light Sources and Probes

ASD offers several accessories for:

- Delivering illumination to the sample
- Collecting reflected or transmitted light from the sample
- Transmitting the collected light to the spectrometer

Some accessories, such as the Contact Probe, perform all of these functions. Other accessories, such as some of the optional foreoptic probes, rely on other illumination sources.

Accurate analysis of collected spectra requires an adequate level of incident light on the sample. The Contact Probe and MugLite accessories are designed for stability and accuracy and provide a controlled source of incident light. They are powered by the accessory power port of the AgriSpec spectrometer. Characteristics of the light source are eliminated when the baseline spectrum is applied to each raw sample spectrum.

Notes:

Chapter 2 Setup

Inspect the shipping container and take careful notes regarding any damage that might have occurred during shipping.

Note: Save all packing materials, foam spacers, and paperwork for possible future use.



Figure 2-1 AgriSpec spectrometer in carrying case.

Prepare a clear space on a sturdy bench or counter, ideally, a space near a wall-current receptacle and the controlling computer.

Carefully open the shipping container following all instructions and orientation labels on the container.

2.1 Unpacking the Instrument

Step 1 Carefully remove the instrument from its shipping case.

Note: During operation, the unit will work correctly regardless of the physical orientation, such as sideways in a specially designed Ergonomic Pro-Pack.



Figure 2-2 AgriSpec spectrometer front-view showing fiber connection and power output.

Step 2 The computer is located directly below the instrument between the foam inserts.

Step 3 Under the computer compartment, there is a circular cut-out. Carefully remove the fiber-optic cable located below the foam insert at the bottom of the case.

The remaining portion of the case is designed to hold the various accessories that may have purchased with your instrument. There are specific locations for the Contact Probe, Hi-Bright Muglite, Spectralon, battery pack, and miscellaneous power cables.

WARNING! The fiber cable should never be stored with a bend of less than a 5" diameter for long periods of time, because the cable can be damaged with undetectable longitudinal fractures that can cause light leakage and weaker signals.

2.2 Power Supply

The AgriSpec spectrometer requires input power to be 12 VDC (60 W). It does not contain an internal power supply to convert AC voltage to DC, nor does it have internal batteries.

The three options for providing the spectrometer with the appropriate DC voltage are:

- An external power supply that converts 100 VAC to 240 VAC (50 Hz or 60 Hz) to 12 VDC (shown in [Figure 2-3](#)).
- An external battery pack that delivers 12 VDC.
- An external adapter for vehicle cigarette lighters (sold separately and should not be used with vehicle engine running).

Caution! *Use only ASD approved power supplies or connectors to power the instrument.*



Figure 2-3 AgriSpec spectrometer power supply and cables.



Figure 2-4 Portable battery pack .

Both the external AC power supply and the DC battery pack have the appropriate power connector (Figure 2-5) to connect to the AgriSpec spectrometer on the back panel (Figure 2-6).



Figure 2-5 AgriSpec spectrometer power connector which plugs into the instrument.

Insert power (battery or power supply) into the receptacle on the back panel labeled “**INPUT 12 VDC 60 WATT**”. Lock the connector in place by twisting the barrel end of the connector clockwise.



Figure 2-6 AgriSpec spectrometer back panel showing Ethernet connection, power switch, and power input jack.

2.3 Ethernet Cables

For Information about configuration of the Ethernet card and cable please refer to the *Installation Guide*. Only use shielded Ethernet cables to reduce electromagnetic interference with the instrument.

Note the location of the computer interface connection. [Figure 2-6](#) shows the back panel containing the Ethernet port, which is labeled “**RJ45/ETHERNET**”.

With the AgriSpec spectrometer and the instrument controller fully powered-down and turned-off, plug the shielded cross-over Ethernet computer cable (as shown in [Figure 2-7](#)) into the Ethernet ports of the AgriSpec spectrometer and the instrument controller.



Figure 2-7 Shielded cross-over Ethernet cable.

2.4 Accessory Power Port

The accessory power port is located on the front of the unit, as shown in [Figure 2-8](#). This port is where external accessories, such as the Contact Probe or Source Probe, attach to receive power from the spectrometer.



Figure 2-8 Enhanced view of the front panel with the fiber optic cable connector, port for remote trigger, and accessory power port for probe.

The accessory power port is provided for optional ASD approved accessories. Attach the large connector end of the cable into the accessory power port of the front panel. Attach the other end of the power cable into the Contact Probe, Hi-Brite MugLite, or other ASD approved accessory.

Please contact ASD for information about these accessories.

WARNING! Never attempt to use unauthorized accessories on the accessory power port



Figure 2-9 Cable used from the AgriSpec spectrometer accessory power port to the Contact Probe, MugLite, or other authorized ASD accessory.

2.5 Remote Trigger

An external trigger is standard with the AgriSpec spectrometer.

The LEDs on the trigger are bright enough to be seen even on a sunny day. The LEDs are lit up when the trigger is pushed (briefly) and turned off when the capture is complete. [Figure 2-10](#) shows the trigger switch with its LEDs.

The trigger plugs into the spectrometer near the fiber optic cable, as shown in [Figure 2-11](#).



Figure 2-10 External trigger.



Figure 2-11 External trigger cable plugged into the AgriSpec spectrometer.

The trigger comes with several pieces of mated velcro, which are intended for use with several accessories. Place a piece of hook velcro onto the trigger; place pieces of mesh velcro onto the accessories.

Note: Accessories should have pieces of mesh velcro placed on both sides so that the trigger can be used right- or left-handed.

Chapter 3 Usage and Maintenance

For personal safety and to prevent damage to the instrument, please review the following safety precautions. All operators should be familiar with this information.

3.1 Light Sources

Use only light sources supplied by ASD Inc. Light sources supplied by ASD are designed to provide levels of illumination and stability of output that complement the performance of ASD's instrumentation.

- Never touch the light source bulb and avoid contact with hot metal components near the bulb! Heat transfer from the light source may make these metal components uncomfortably warm to the touch.
- Prevent dirt and oils from contacting the bulb and reflector.
- Do not touch the glass envelope of the bulb. Oils on the hand may transfer to the glass envelope, which can significantly affect the bulb temperature and corrupt important operating physics of the bulb.

Note: Please refer to the *ASD Accessories Manual* or the instructions for that specific accessory for detailed information.

3.2 Access and Ventilation

When used inside, provide adequate room ventilation for the spectrometer. Insufficient ventilation can result in overheating of the spectrometer and subsequently corrupted data and possibly physical damage to the spectrometer.



Do not cover the vents of the instrument.

Do not place objects on the unit or its power supplies. Prevent objects from obstructing ventilation slits.

Keep objects and spills from entering or falling onto the instrument, power supplies, and software disks.

3.3 Cleaning

- Step 1 Turn off the unit.
- Step 2 Disconnect from all power.
- Step 3 Allow the unit to cool down.
- Step 4 Clean unit with a slightly damp cloth and mild soap.

Note: Be sure all soap residue is removed and all surfaces are dry before use.

3.4 Annual Maintenance

ASD recommends that the instrument be serviced once a year. This will ensure the proper function of the instrument. Annual maintenance is covered under ASD warranty or the extended service contract. Annual maintenance can be purchased separately. If you are interested in purchasing annual maintenance or an extended service contract, please contact your sales representative.

3.5 Returning Instrument to ASD for Service

In order to return the instrument to ASD for maintenance or repair, a Return Merchandise Authorization (RMA) must be issued by ASD technical support. The RMA includes scheduling details, contact information, shipping instructions, as well as a brief description of the maintenance or repair requirements. Please contact the technical support department at 303-444-6522 x-144 or support@asdi.com.

Chapter 4 *Fiber Optic Interface*

ASD offers a variety of fiber optic cable lengths to interface with the AgriSpec spectrometer. These fiber optic cables are used for reflectance and transmittance measurements.

Fiber optic cables should be loosely coiled and stored in the compartments provided. All accessories should be stored in the designated compartments of the carrier or in a separate protective case.

4.1 *Fiber Optics Cables*

Two fiber optic cables are included with the spectrometer: an external cable and an internal cable (built into the instrument).

- The built-in fiber optic cable is made up of fifty-seven (57) randomly distributed, ultra low-OH, silica glass fibers. Nineteen (19) of these fibers are 100 micron and are distributed to the VNIR region. The remaining thirty-eight (38) fibers are 200 micron and are evenly divided between the two SWIR regions.
- The external fiber optic cable is made up of forty-four (44) randomly distributed, ultra low-OH, silica glass fibers that are 200 micron. This cable has a Numerical Aperture (NA) of 0.22.

The external fiber optic cable length is typically 1 meter. Additional external fiber optic cables can be ordered in a variety of lengths. However, as the length increases, the cable's transmission decreases at wavelengths below 400 nm and above 2000 nm.

External fiber optic cables should be coiled and stored in the compartment provided in the shipping case. All accessories should be stored in the designated compartment of the case or in a separate protective case.

The external fiber optic cable are protected with a metal spiral inside the black PVC cable casing. If there are kinks in the cable, the fibers are not necessarily damaged. If, however, the cable has been damaged so severely that the protective metal spiral can be seen, the chances are high that the fibers have been damaged. Each broken fiber results in a ~5% loss of response.

WARNING! The fiber optic cable should never be stored with a bend of less than a 5" diameter for long periods of time.

The fibers can be damaged by coiling the cable up too tightly. If left in a tight coil for longer than a week, the fibers are likely to develop longitudinal fractures that will not be detectable. These fractures in the fiber will cause light leakage, resulting in a weaker signal. The fiber optic cables should be stored by placing them loosely within an appropriate compartment of the carrying case or in their own protective case.

Tips on care for the fiber optic cable:

- Do not pull or hang the spectrometer by the fiber optic cable.
- Do not use wires, ties, or clamps to tightly attach the fiber optic cable to objects, because these might pinch or penetrate the protective jacket thereby damage the fibers inside.
- Avoid whipping the fiber optic cable, dropping it, or slamming it into objects, because this can cause fractures to the glass fibers.
- Avoid twisting the fiber optic cable, because such forces may cause fractures to fibers.

While the tip of the fiber optic cable is not particularly susceptible to damage, a tip cover is recommended to protect against abrasion and exposure to contamination. Replacement covers can be made by cutting pieces of eighth-inch shrink tubing to about 1.5" lengths and shrinking them onto the fiber cable tip. The covers will slide on and off the cable easily.

4.2 Fiber Optic Cable Interface

All fiber optic cables come with a standard SMA style 905 connector, as shown in [Figure 4-1](#)

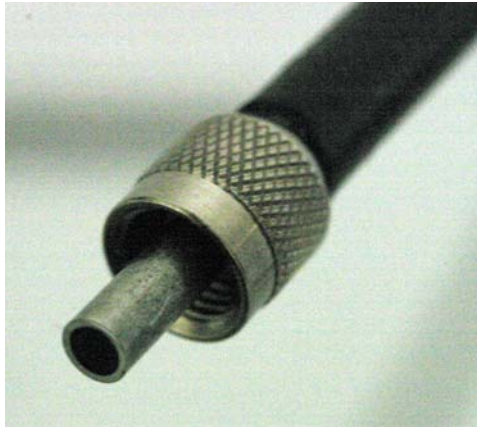


Figure 4-1 Connector for the fiberoptic cable.

In order to achieve the best response from ASD instruments it may be necessary to adjust the fiber optic interface between the external fiber optic cable (shown in [Figure 4-2](#) and [Figure 4-3](#)) and the instrument connection.



Figure 4-2 View of the fiberoptic bundle matcher and the fiberoptic port.

Note: The fiber optic bundle matcher has a standard 1.7 mm fiber optic bundle input.



Figure 4-3 View of the fiberoptic cable, fiber bundle matcher, and fiber connection. Screw the fiber bundle matcher into the fiber connection as far as possible.

ASD external fiber optic cables most commonly come in one of the configurations listed in Figure 4-4. The millimeter size (mm) is the internal diameter of the fiber optic bundle.

The AgriSpec spectrometer is almost always used with the 1.7 mm external fiber optic cable diameter.

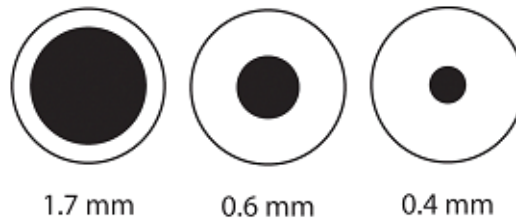


Figure 4-4 Examples of the external fiberoptic cable diameters.

4.3 Matching Fiber Bundles

Now that there is an understanding of how the external fiber optic cables and the fixed fiber optic bundle matcher are configured, they need to be combined. The three scenarios in Figure 4-5 show how the fiber optic bundles are matched.

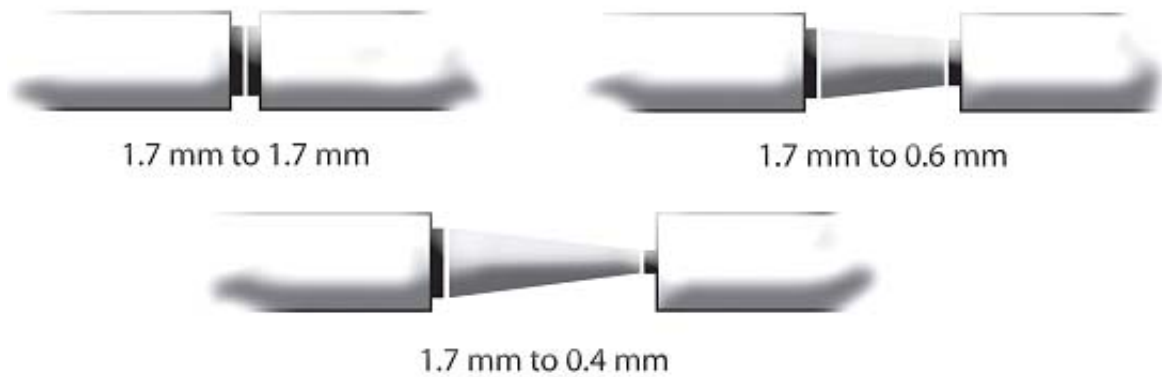


Figure 4-5 Matching fiber bundles to the AgriSpec spectrometer fiber port.

Note: The optimal setting for this fiber optic bundle matcher is adjusted at the ASD factory. If the customer has ordered several fiber optic configurations, it may be necessary to adjust this bundle matcher in order to achieve the highest signal to noise ratio and best instrument response. Changing the bundle matcher can effect the performance of the spectrometer.

In order to match the bundles:

- 1 Unlock the hex-nut bolt at the front of the bundle matcher on the spectrometer with the 3/8 in hex-nut driver (included with the instrument.)
- 2 The AgriSpec spectrometer should be set up with the adjustment barrel of the matcher fully threaded into the body of the matcher any time a full 1.7 mm bundle is used (which is almost always).
- 3 With the ASD software application (RS³ or Indico) running and the fiber optic cables attached, hand screw the fiber optic bundle matcher until the optimal signal is achieved. *Do not over twist the fiber optic bundle!*

The optimal signal will be achieved by watching the RAW DN (DN = digital numbers) in the ASD application as the fiber optic bundle matcher is being hand adjusted.

Note: The AgriSpec spectrometer should be set up with the adjustment barrel of the matcher fully threaded into the body of the matcher any time a full 1.7 mm bundle is used.

- 4 Once the optimal signal is achieved, lock down the fiber optic bundle matcher. Tighten the 3/8 hex-nut by turning the wing nut against the face of the instrument. Use the hex wrench to hold the 3/8 hex-nut in place while you tighten the wing nut.

When the bundle matcher is removed, clean the fiber optic cable and the exposed fiber optic on the spectrometer. This can be done by using de-ionized water, cleaning the end, and then gently wiping the excess off. Canned air to blow off any dust that may have accumulated can be used.

4.4 LED Check for Fiber Optic Cable

The AgriSpec spectrometer allows you to perform a visual verification of the internal fiber optic cable using built-in LEDs, a magnifier placed on the fiber optic bundle matcher, and the **Fiber Check** software.

WARNING! The **Fiber Check** utility produces rapid flashing lights in the SWIR 1 and 2 region's optical fibers. If you are susceptible to epileptic seizures, exercise caution or avoid using the **Fiber Check** utility.

- Step 1 Remove any external fiber optic cable from the spectrometer.
- Step 2 Remove the bundle matcher. Remove the bundle matcher. Use the 3/8 inch hex-nut driver to loosen the bundle matcher if necessary.
- Step 3 Remove the Fiber Optic Port using the 9/16 nut driver (included with the instrument.)

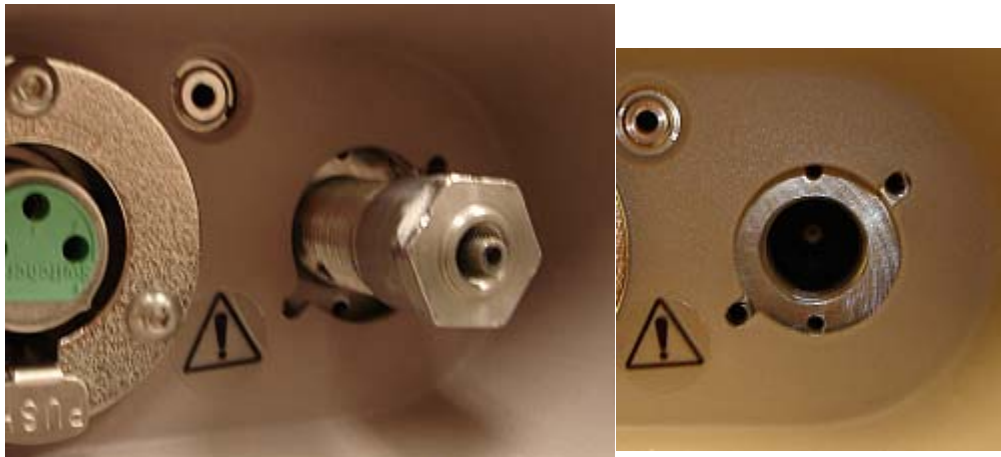


Figure 4-6 Fiber Optic Port

- Step 4 Carefully insert the tip of the magnifier until it butts up against the internal fiber optic, as shown in [Figure 4-7](#).

Note: Be careful not to apply any load or torque to the magnifier which could break the tip off.

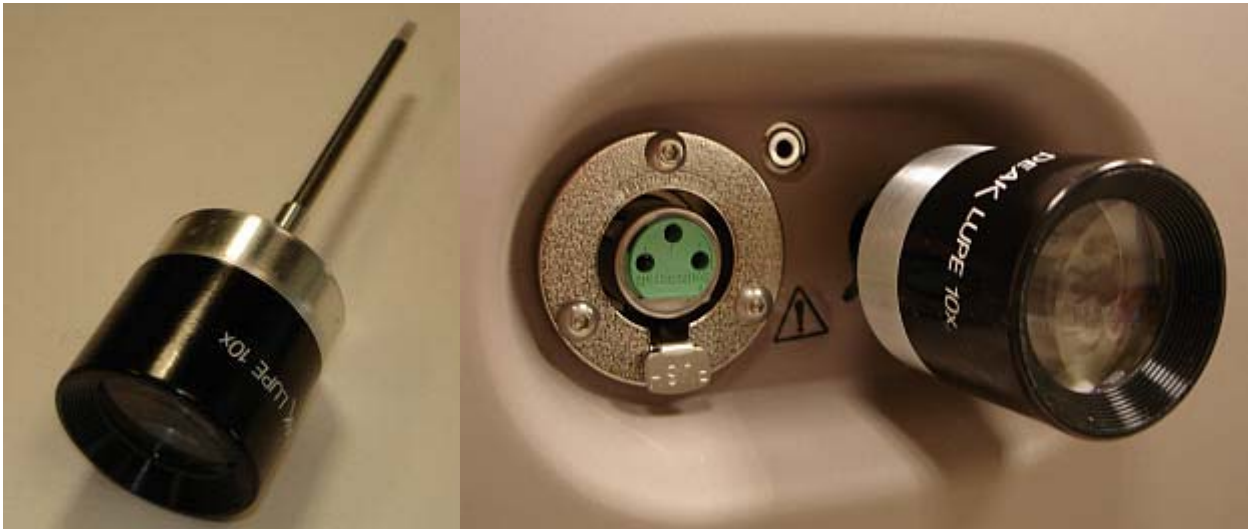


Figure 4-7 Magnifier for testing the internal fiber optic cable.

- Step 5 On the instrument controller, exit any ASD applications that might be running and communicating with the unit. Such applications include RS³ and Indico.
- Step 6 Ensure that the spectrometer is turned on.
- Step 7 Start the **Fiber Check** application from the **Start** menu under **All Programs->ASD Programs->Indico Pro Tools->Fiber Check**.

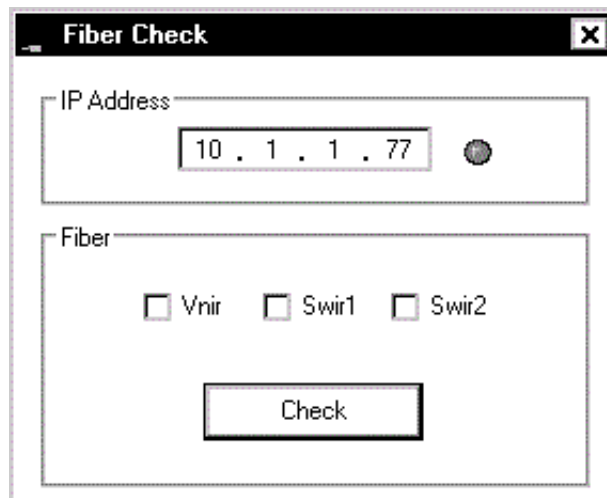


Figure 4-8 **Fiber Check** Program

- Step 8 Ensure that the IP address configured is the one for the spectrometer.

Step 9 Use the mouse to select the checkbox for the LED to turn on: VNIR, SWIR1, or SWIR2. One or more may be selected at once.

Note: If a range is selected that the spectrometer does not have, that particular LED will not be available to turn on. Nothing will be harmed.

Step 10 Select the **Check** button which turns on the selected LEDs.

Step 11 Look through the magnifier to see which fibers light up.

- Count the number of fibers that show light. Refer to [Details about the Fiber Optic Cable and LEDs](#).
- If applicable, use the checkboxes and the **Check** button shown in [Figure 4-8](#) to turn on and off different LEDs to help determine which range might be affected.

Note: The RED LED for VNIR may be hard to see when the other ranges are enabled.



Figure 4-9 Internal spectrometer LED shining to test internal fiber optic cable.

Step 12 When you are finished, shutdown the **Fiber Check** application.

Step 13 Carefully remove the magnifier and re-attach the bundle matcher and fiber optic port using the 9/16 nut driver.

Step 14 Inspect the external fiber optic cable disconnected from the instrument.

- 1 Unscrew the lens from the magnifier and place it onto the short magnifier assembly



Figure 4-10 Short and Long Magnifier Assemblies

- 2 Attach the magnifier to one end of the fiber optic cable.
- 3 Point the other end of the fiber optic cable close to a light.
- 4 Observe which fibers transmit light.

Details about the Fiber Optic Cable and LEDs

- The LEDs will only turn on for the ranges installed in the spectrometer. Selecting one that the spectrometer does not have will not cause a problem.
- Each range has a different LED color so that you can see the fibers in the cable that are associated with it: RED for VNIR; WHITE for SWIR1; GREEN for SWIR2.
- The fiber bundle of the internal cable for each range contains 19 fibers.
- The fiber bundle of the external cable contains 44 fibers.
- Each broken *internal* fiber results in an approximate 5% loss of response in that particular range (VNIR, SWIR1, or SWIR2). Each broken *external* fiber results in a few percentage loss of response over the entire range. It is hard to quantify how a broken fiber in the external cable affects the response of any given range, each time the external cable is connected to the spectrometer, the alignment with the fibers of the internal cable will most likely be different.


The instrument can be successfully used with a few broken fibers in each range, although with a reduction in signal strength.

For severely damaged internal fiber optic cables, send the instrument back to ASD Inc. in its carrying case for repairs. For severely damaged external fiber optic cables, order replacements from ASD Inc.

Notes:

Chapter 5 Batteries

The specifications of the battery are as follows:

Type:	Gel Cell Electrolyte
Rating:	12 V 9 amp hour
Life:	Over four (4) hours using a full-range AgriSpec spectrometer with a Contact Probe in an ambient environment. Nine (9) hours using a VNIR-only AgriSpec spectrometer with accessory not powered by the instrument in an ambient environment.
	Recycle batteries and do not dispose of as general waste.

To run the instrument with battery power, plug the battery cable from the battery to the instrument power connector (back of instrument).

Note: How long the spectrometer will run with a fully charged battery depends on many variables including battery age, instrument configuration, environment, and accessories powered by the accessory power port.

5.1 Battery Care

Proper care of the Gel Cell Electrolyte rechargeable batteries ensures a long life.

The batteries are rated to give 200 to 300 uses when they are fully drained (discharged to 9 Volts) each use. Letting the battery discharge under 9 Volts will shorten the battery life. If the battery is only used for 50% rated capacity up to 500 uses are possible. In the same way using only 30% capacity can result in up to 1,000 uses.

Note: There is no such thing as memory effect on these cells. Only use them for the time needed. When finished with the batteries, recharge them fully and then store them away until next use.

Recharge a fully discharged battery as soon as possible (usually within 24 hours). Uncharged batteries will be incapable of accepting a charge if the period of time between charges is too long. A partially discharged battery does not need charging immediately. However, if the battery is to be stored away, first charge it to full capacity. When 100% charged, these batteries can be stored up to six months, at that time perform a topping charge before use. Always store the battery in a cool, dry location.

WARNING! Do not use batteries other than those supplied by ASD. Do not use ASD batteries in a manner unauthorized by ASD. Using improper batteries or improper use of ASD batteries could result in bodily injury or damage to the instrument.

5.2 Battery Charger

The battery charger specifications:

Type:	External Desktop
AC Input:	90-240 VAC, 50/60 Hz
DC Input:	12 V 1.5 A
Charge Time:	under 8 hours for a fully discharged 9 AH battery

Review the charger instruction manual for details.

LED Status indicator:

- The LED will be ORANGE during charge.
- The LED will turn GREEN when charge is completed.

The battery charger is designed for indoor use and should not come in contact with water or dust.

Note: Do not unplug the charger and leave the batteries in the charger. Remove the batteries first, or they will discharge through the charger.

Constant overcharging or undercharging will shorten battery life.



The AgriSpec spectrometer will stop collecting spectra when the battery drains to about 9.7 volts and the software will post a “Spectrometer Isn't Ready” message. However, the instrument and attached probe will continue to draw current if power is left on. Power down the instrument and remove probe as soon as possible when the above message occurs.

5.3 Application Battery Indicator

The RS³ and Indico applications from ASD have a **Power Status**. It is located at the bottom of the main window next to the **Connection Status**. The voltage level of the spectrometer will be reported in the balloon type window.

- 11-12 Volts is considered good (full).
- <11 Volts issues a warning.

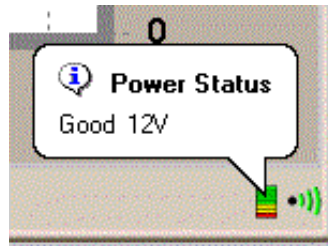


Figure 5-1 Example of a **Power Status**

Notes:

Chapter 6 *Field Measurements*

Spectrometry is the quantitative measurement of reflectance, transmission, or absorption. It involves the collection of accurate spectra and requires an awareness of the influences of:

Accurate and repeatable material analysis or identification requires the accurate and repeatable measurement of reflectance, transmission, or absorption. It requires an awareness of the influences of:

- Sources of illumination.
- Instrument field-of-view.
- Sample viewing and illumination geometry.
- Instrument scanning time.
- Spatial and temporal variability of the sample characteristics.

Many of these parameters are controlled when using one of the ASD standard sampling interfaces (e.g., MugLite or Contact Probe).

Spectral measurements are desirable for many reasons.

- Spectra of target materials allows for more precise image analysis and interpretation.
- Spectroscopy can perform feasibility studies to understand if and how a process or material of interest can be detected using remote sensing.
- Field spectra can be used to make direct material identifications in the field rather than collecting samples for later laboratory analysis.

The first step in the development of a field experiment is the definition of the overall experimental design. Unfortunately, the formulation of an appropriate experimental design is not always obvious.

Issues such as the timing of the data collection, spatial scale of the field measurement, target viewing and illumination geometry, and the collection of ancillary data sets must be considered in light of the objectives of the study. The lack of appropriate ancillary data sets often makes previously collected data sets unusable for a new application.

Frequently, the experimental design must be modified to account for the characteristics of the available instrumentation.

6.1 Illumination

In order to determine the reflectance or transmittance of a material, two measurements are required:

- The spectral response of a reference sample.
- The spectral response of the target material.

The reflectance or transmittance spectrum is then computed by dividing the spectral response of the target material by that of a reference sample.

Using this method, all parameters which are multiplicative in nature and present in both the spectral response of a reference sample and the target material, are ratio-ed out, such as:

- The spectral irradiance of the illumination source.
- The optical throughput of the field spectrometer.

Note: An inherent assumption when determining the reflectance or transmittance of a material in the field is that the characteristics of the illumination are the same for the reference and target materials. *Variability of the illumination characteristics* between the time the reference and target materials are measured *will result in errors in the resultant spectra.*

6.2 Vegetation

The absorption features seen in vegetation spectra are all related to organic compounds common to the majority of plant species.

Thus, the information about a plant canopy is contained in the relative intensity of the various absorption features rather than in the presence or absence of a specific absorption feature.

The major spectral absorption features can be attributed to:

- water
- plant pigments
 - » chlorophyll
 - » zanthophyll
 - » carotenoids

Other, minor, absorption features are attributable to other chemical components; these include:

- cellulose
- lignin
- proteins

- starches
- sugars

Non-photosynthetic components of the canopy have spectra which are dominated by absorption features attributed to lignin and cellulose.

6.3 Rocks, Soils, and Man-Made Materials

The shape of the spectral signature of rocks and soil tends to be invariant with respect to the viewing geometry, unlike vegetation. The overall brightness of the observed spectrum does change with illumination and viewing geometry due to changes in the amount of shadow in the field-of-view of the spectrometer.

Absorption features in the spectra of rocks and minerals are due to the presence of specific molecular groups and are often diagnostic of the minerals present in the sample.

6.4 Suggested Set-up

6.4.1 RS3 Application

- Step 1 Follow the instructions for plugging in the instrument and starting the RS³ software.
- Step 2 Attach an appropriate accessory. Remember that the AgriSpec spectrometer requires an external (artificial) light source.
- Step 3 Open the pull down menu **Control -> Adjust Configuration [Alt+C,C]** in the RS³ application.
- Step 4 Set the white reference number of scans to 30, and dark current averages to 60 scans.
- Step 5 Set the spectrum average (or sample count) to 30 for indoor or artificial illumination.
- Note: The actual spectrum average will be determined by striking a compromise between noise reduction through averaging the spectra and the time desired for each spectrum collection. For instance, if you are using the instrument in the field, are walking a large area, and are making frequent spectral readings, you will want a shorter average setting than if you are collecting spectra in-situ and desire the cleanest spectra possible.
- Step 6 **Optimize** the instrument to the Spectralon.

Spectralon is an example of a reference standard to take a white reference against. (Labsphere, Inc., Shaker St., PO Box 70, North Sutton, NH 03260-0070, PH: 603-927-4266, FX: 603-927-4694, E-mail: sales@labsphere.com, Website: www.labsphere.com). Spectralon is also available from ASD.

- Step 7 Take a **White Reference** (or baseline reference).
- Step 8 Observe samples.

6.4.2 Indico Application

- Step 1 Follow the instructions for plugging in the instrument and starting the Indico software.
- Step 2 Attach an appropriate accessory. Remember that the AgriSpec spectrometer requires an external (artificial) light source.
- Step 3 Open a new project either by selecting **File->New Project** from the pull-down menu.
- Step 4 Type in a name for this project and press the **OK** button.
- Step 5 Select **Spectrum->Sample Count/Average** from the pull-down menu.
- Step 6 Set the **Instrument Sample Count** to 30 for indoor or artificial illumination.

Note: The actual spectrum average will be determined by striking a compromise between noise reduction through averaging the spectra and the time desired for each spectrum collection. For instance, if you are using the instrument in the field, are walking a large area, and are making frequent spectral readings, you will want a shorter average setting than if you are collecting spectra in-situ and desire the cleanest spectra possible.

The project screen will return, and the sample count will be displayed in the status box at the bottom of the graph.

- Step 7 Ensure that the light source is on and the probe input end is pointed at the white reference panel.
- Step 8 Select **Spectrum->Take a Baseline [Alt+S, B] [Ctrl+R]**. When prompted, click **Yes, take a new white reference**.
 - The AgriSpec spectrometer optimizes the detector sensitivities for the probe and light source currently being used.
 - The dark offset and white reference will also be measured and saved.
 - Status bars will indicate each process.

When optimization and white reference are complete, you should see a straight baseline across the Project Graph at 1.00 (100%).

- Step 9 Select the **Display->Reflectance [Alt+D, R]** for the reflectance mode.

- Step 10 Place the sample on the white reference panel and the spacer end in contact against the sample.
- This assumes that your reflectance probe has appropriate spacing or spacers on the end.
- Step 11 Press the Space Bar once or select the **Spectrum->Take a Scan [Alt+S, T]**.
- When the AgriSpec spectrometer finishes recording and averaging the Spectra, the graph in the project window will be updated to show the reflectance spectrum of the sample.
- Step 12 Observe samples.

6.5 White Reference

A material with 95-99% reflectance across the entire spectrum is called a *white reference panel* or *white reference standard*.

The raw measurement made by the spectrometer is influenced by both the sample and the light source. An independent measure of the light source illumination on a reference of known reflectance is required to calculate the reflectance of the sample. The use of a white reference standard with near 100% reflectance simplifies this calculation.

The ASD application software, such as RS³ and Indico, can calculate the ratios for reflectance or transmittance of the material being sampled by the AgriSpec spectrometer using the white reference as the standard.

Spectralon from Labsphere is the white reference standard that is very suitable for the VNIR and SWIR spectral ranges of ASD instruments.

Spectralon is made of polytetrafluoroethylene (PTFE) and cindered halon. It has the characteristic of being nearly 100% reflective within the wavelength range of 350 nm to 2500 nm. A Spectralon white reference scatters light uniformly in all directions within that wavelength range.

6.5.1 Spectralon Reflectance Data

Table 6-1 shows the published reflectance data for an uncalibrated SRM-990. When SRM-990 is used as the white reference for a reflectance measurement with the AgriSpec spectrometer, an even closer reflectance value for the sample can be calculated.

Table 6-1 Labsphere's published reflectance data for uncalibrated Spectralon (+/- 0.5%)

Wavelength (nm)	SRM-990 "Uncalibrated" Spectralon Reflectance
250	0.950
300	0.985
400	0.990
500	0.991
600	0.992
700	0.992
800	0.991
900	0.991
1000	0.993
1100	0.993
1200	0.992
1300	0.992
1400	0.991
1500	0.991
1600	0.991
1700	0.988
1800	0.989
1900	0.981
2000	0.976
2100	0.953
2200	0.973
2300	0.972
2400	0.955
2500	0.950

Reflectance = Energy reflected from target/energy incident on target

6.5.2 Maintaining Spectralon References

Spectralon is an optical standard and should be handled in much the same way as other optical standards. Although the material is very durable, care should be taken to prevent contaminants such as finger oils from contacting the material's surface. Always wear clean gloves when handling Spectralon.

To clean a lightly soiled Spectralon white reference

If the material is lightly soiled, it may be air brushed with a jet of clean dry air or nitrogen.

WARNING! DO NOT use Freon.

To clean a heavily soiled Spectralon white reference

Sand the Spectralon material under running water with a 220-240 grit waterproof emery cloth or sandpaper until the surface is totally hydrophobic (water beads and runs off immediately).

- Step 1 Use a flat surface, such as a thick, flat piece of glass.
- Step 2 Place the glass into the sink.
- Step 3 Place 220 grade wet sandpaper onto the glass.
- Step 4 Gently move the Spectralon reference in a figure 8 motion on the sandpaper, using water as needed to wash away the thin layer that is sanded off.
- Step 5 Blow dry with clean air or nitrogen or allow the material to air dry.
- Step 6 If the material requires high resistance to deep UV radiation, the piece should be prepared as above, then either of the following two treatments performed.
 - 1 Flush the Spectralon piece with >18 milli-ohm distilled, deionized water for 24 hours.
 - 2 Vacuum bake the Spectralon piece at 75° C for a 12 hour period at a vacuum of 1 Torr or less. Then purge the vacuum oven with clean dry air or nitrogen.

WARNING! Do not use oils or soaps to clean the Spectralon white reference.

6.5.3 White Reference Procedures

Baseline references, or white references, should be collected approximately every 15 minutes while the instrument is warming up; thereafter, every hour or so is sufficient. The Spectralon puck should be used when optimizing and taking a white reference measurement.

The AgriSpec spectrometer should be re-optimized for:

- Light changes.
- Temperature changes.
- Whenever accessory probes are changed.

Note: Environmental conditions can change rapidly or slowly. It all depends on clouds, wind (affecting temperature), instrument warm up time, etc.

Appendix A Troubleshooting

A.1 Common Communication Fixes

To fix many communication errors (particularly if the instrument has been functioning at some point), power cycle the instrument and/or the instrument controller.

The sequence to use will vary depending on the computer manufacturer. Either:

- Leave the computer on. Turn off the instrument. Wait for 10 seconds. Turn the instrument back on.
- Or, turn off the computer and the instrument. Turn on the instrument. Then turn on the instrument controller.

For the best performance, you should ensure a clear line-of-sight between the spectrometer and instrument controller. Obstructions or radio frequency interference in close proximity can cause loss of communication or significantly shorten the communication range.

The AgriSpec spectrometer uses industry standard components. You should see the same general connection speed and distance capabilities as other wireless devices.

A.2 Does Not Connect to the Spectrometer

- Refer to section [A.1, Common Communication Fixes](#).
- Check that the Ethernet cable is securely inserted into the spectrometer and host computer. Check that the Ethernet LED is on.
- For an isolated network, check that the Ethernet cable is a cross-over Ethernet cable.
- For an established network, check that the Ethernet cable is a standard Ethernet cable.
- Check that the IP Address is in the same range or subnet as the ASD spectrometer. The same subnet means that the first three octets of the IP address (xxx.xxx.xxx.___) match the spectrometer and the computer.

- Do a ping test to make sure the spectrometer is responding.
 - » Open up a command window by selecting **Start->Run**
 - » Type **cmd** in the **Run** window.
 - » Select **OK** to open the window.
 - » For an Ethernet connection, type: ping 10.1.1.11
 - » For a wireless connection, type: ping 10.1.1.77

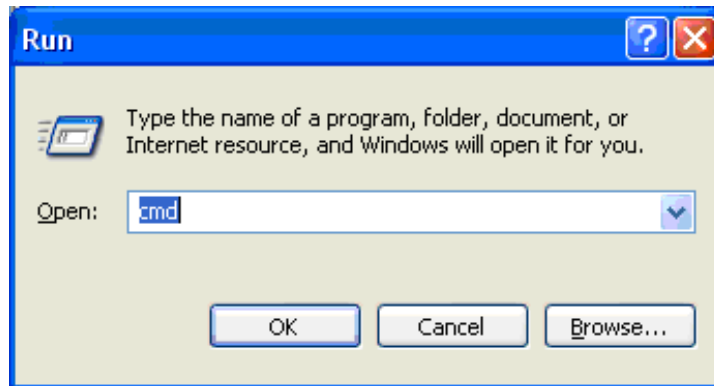


Figure A-1 Run window with **cmd**.

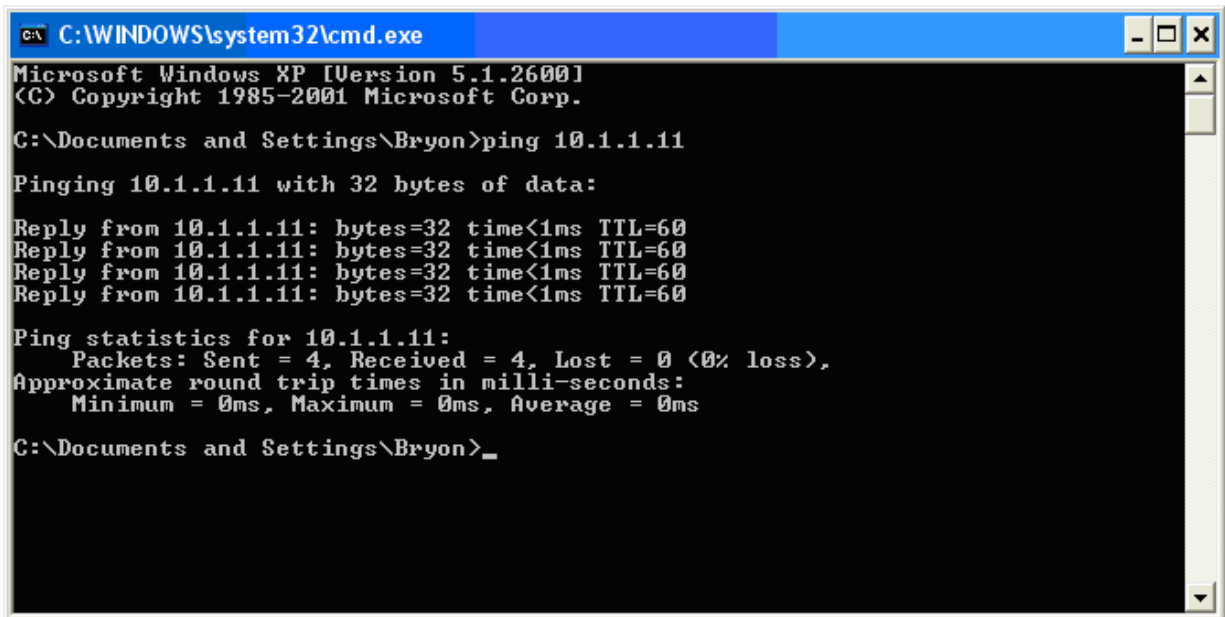


Figure A-2 Successful result of a **ping** operation.

The default IP Address for the ASD spectrometer is:

- 10.1.1.11 for the Ethernet interface, or
- 10.1.1.77 for the wireless interface.

The default subnet mask is 255.255.255.0.

The computer's Ethernet adapter or wireless adapter must have a unique IP address in the same range as the spectrometer, such as 10.1.1.x where **x** is a unique number. The subnet mask must also be the same as the spectrometer, such as 255.255.255.0.

A.3 Does Not Connect Wireless

- Refer to section [A.1, Common Communication Fixes](#).
- Check to ensure that the wireless adapter is in the same range as the ASD spectrometer.

The default IP Address for the ASD spectrometer is:

- » 10.1.1.11 for the Ethernet interface, or
- » 10.1.1.77 for the wireless interface.

The default subnet mask is 255.255.255.0.

The computer Ethernet adapter or wireless adapter must have a unique IP address in the same range as the spectrometer, such as 10.1.1.x where **x** is a unique number. The subnet mask must also be the same as the spectrometer, such as 255.255.255.0.

- Check to ensure that the wireless adapter is connected to the spectrometer. (For Windows XP)
 - » Select **Start->Network Connections**.
 - » Right click on the wireless adapter.
 - » Select **View Available Wireless Networks**.
 - » Ensure that the ASD spectrometer shows up in the list of wireless networks. If the spectrometer is not connected,
 - » Select the spectrometer.
 - » Select the **Connect** button.
 - » If you are using encryption, then make sure you are using the same WEP key in the wireless adapter as the ASD spectrometer.
 - » Power cycle the spectrometer.

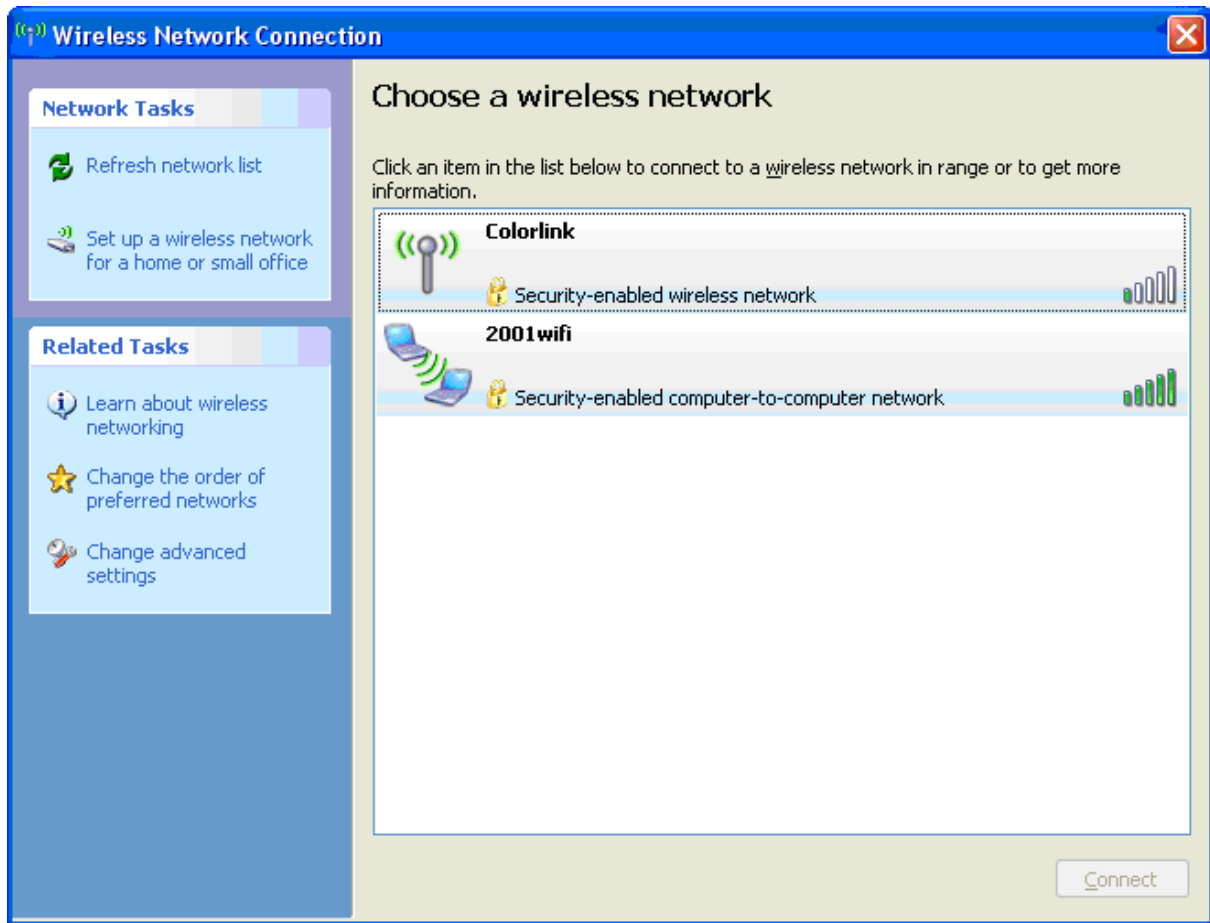


Figure A-3 Wireless Network Connection.

A.4 Does Not Connect to the Access Point

- Refer to section [A.1, Common Communication Fixes](#).
- Check to ensure that the wireless interface on the spectrometer is not set to **ad hoc**.
- Check to ensure that the **SSID** of the spectrometer is the same as the access point.
- If you are using encryption, then ensure that you are using the same **WEP** key in the spectrometer as the access point.
- Move the access point, spectrometer, and instrument controller (computer) to the same room to test the connection.

A.5 Spectrometer Loses its Wireless Connection

- Refer to section [A.1, Common Communication Fixes](#).
- Try to minimize the RF noise in your environment. RF noise interference can come from: 2.4 Ghz cordless phones, microwaves, monitors, electric motors, ceiling fan, lights, security systems, etc.
- Positioning of the spectrometer and instrument controller (computer) can affect the wireless range. Minimize obstructions between the spectrometer and the instrument controller. Wall, ceilings, doors, buildings, hills, etc. can degrade the signal of the spectrometer and computer.

A.6 Wireless Interface is Disabled in the Netcfg Application

- Refer to section [A.1, Common Communication Fixes](#).
- Check to ensure that the wireless adapter is in the same range as the spectrometer.

The default IP Address for the ASD spectrometer is:

- » 10.1.1.11 for the Ethernet interface, or
- » 10.1.1.77 for the wireless interface.

The default subnet mask is 255.255.255.0.

The computer Ethernet adapter or wireless adapter must have a unique IP address in the same range as the spectrometer, such as 10.1.1.x where **x** is a unique number. The subnet mask must also be the same as the spectrometer, such as 255.255.255.0.

- Unplug the Ethernet cable from the spectrometer.

Notes:

Appendix B Reference Information

B.1 Frequently Asked Questions (FAQ)

B.1.1 What is a spectrometer?

Spectrograph - An optical instrument for forming the spectrum of a light source and recording it on a film. The dispersing medium may be a prism or a diffraction grating. This term was common prior to the digital age. ASD instruments do not use film.

Spectrometer - An optical instrument which uses detectors other than photographic film to measure the distribution of radiation in a particular wavelength region. All ASD instruments are spectrometers.

Spectroradiometer - An optical instrument for measuring the radiant energy (radiance or irradiance) from a source at each wavelength throughout the spectrum. A spectroradiometer is a special kind of spectrometer.

Note: Only the FieldSpec spectrometer is configured as a spectroradiometer. The reason is that the fiber optic cable significantly affects the calibration required to measure radiance. The FieldSpec instrument has a fixed fiber optic cable routed directly through the front panel to the internal spectrometers, while the other ASD spectrometers are designed to detach and even interchange fiber optic cables. To be meaningful, radiance measurements with non-FieldSpec instruments would have to be performed using the **same fiber optic cable** and the **same connector orientation** of the plugged in fiber optic cable **as the factory calibration**.

The SWIR component of the ASD spectrometer is a *scanning spectrometer*, while the VNIR component is an *array spectrometer*.

B.1.2 What are the differences between the ASD spectrometers?

FieldSpec - is a general-purpose spectrometer specifically designed for field environmental remote sensing. The FieldSpec is a spectroradiometer in that it is calibrated to be able to perform radiance and irradiance measurements, as well as reflectance and transmittance. The fiber optic cable is routed through the front panel and is fixed to the internal spectrometers, which provides a superior signal-to-noise ratio than the other ASD spectrometers with detachable fiber optic cables. (The FieldSpec spectrometer can also be specially ordered as a high-resolution instrument, the main TerraSpec feature.)

TerraSpec - is a high-resolution spectrometer specially designed for mining and geology applications. It is ideally suited for performing reflectance spectral measurements of rocks, minerals, and core samples. Its higher resolution is optimized for samples that have high absorption features. The Hi-Brite MugLite and Hi-Brite Contact Probe are designed to work with this spectrometer.

AgriSpec - is a general-purpose spectrometer designed for agriculture and the field analysis of vegetation and soils. Its resolution is ideally suited for performing reflectance and transmittance spectral measurements of samples that have broad absorptions features. This spectrometer can use nearly all of the accessories containing a light source, such as the plant probe and leaf clip.

LabSpec - is a general-purpose spectrometer designed not only for laboratory analysis, but also for the plant floor, receiving, and even field work. It is meant to be portable for multiple uses requiring visible and near infrared reflectance, transmission, or absorbance to analyze and identify materials. The LabSpec is similar to the AgriSpec. The LabSpec 2500 has a 10 nm resolution, while the LabSpec 2600 has a 6 nm resolution.

QualitySpec - is a general-purpose spectrometer specifically designed for material analysis and identification in the laboratory. The QualitySpec features are equivalent to the AgriSpec except that it is not field-portable.

B.1.3 What is remote sensing?

Remote sensing is defined as the art and science of obtaining information about an object without being in direct physical contact with the object.

The ASD spectrometers achieve this using visible near-infrared (VNIR) and short-wave infrared (SWIR) spectra.

B.1.4 How often do I need to optimize?

Optimization has no effect on the data unless the measurement has drifted towards saturation, in which case optimization is mandatory. The RS³ application provides a saturation warning.

B.1.5 How often do I need a (baseline) White Reference?

Inside Use or Constant Lighting Conditions

When using the instrument inside under constant lighting conditions or when using an accessory with its own light source, collect a (baseline) white reference every 10 to 15 minutes while the instrument is warming up and then every hour or so thereafter.

Outside Use

When using the instrument outside, you should collect a new white reference frequently for:

- temperature changes
- accessory changes (like switching probes)

It is important that the light intensity is the same for when the white reference is collected as for the measurements of the sample.

B.1.6 What spectrum average (or sample count) should I use?

When used with a MugLite, Contact Probe, or other accessory, it is usually sufficient to use 10-20 scans in the spectrum averaging (or sample count) for the sample and 20-40 scans for the white reference. The white reference isn't taken as often, which is why you would want to include more scans in its averaging.

The signal-to-noise measurement increases with the square root of the number of scans used in the averaging.

Note: The actual spectrum average will be a compromise between noise reduction through spectra averaging and the time required for each spectra collection. For instance, if you are using the instrument in the field to measure a large number of samples, you will want a smaller number of spectra in the average to decrease the collection time required. If you are collecting spectra in the lab, you'll want to increase the number of spectra in the averaging to obtain the cleanest spectra possible.

B.1.7 What if I want to keep the same settings all day?

Even if you could keep the same settings for the integration time, gain value, and offset value for the entire day, you probably wouldn't want to.

Most of the measurements made by the instrument are ratios where such values get cancelled out and become irrelevant.

It is far more important to keep the instrument from saturation and to ensure the (baseline) white reference is fresh.

B.1.8 What are the units of radiance?

Radiance is the radiant power per unit source area per unit solid angle. This is measured in Watts per square meter per steradian ($\text{W}/\text{m}^2/\text{sr}$).

The AgriSpec spectrometer does not measure radiance, because this requires radiometric calibration files which are dependent on the fiber optic cable and its connection to the instrument. (The AgriSpec spectrometer can be used with different accessories and fiber optic cables.)

The AgriSpec spectrometer measures only reflectance and transmission measurements.

B.1.9 Can I post-process my data?

Yes. ViewSpec is one of many applications that can post-process your data. Various software packages are available for chemometric model. The spectral data can be imported into many different applications.

The complete specification of the ASD file format is available upon request.

B.1.10 Why do I see oscillations (sine wave) in my data?

Your light source may use AC power. A single SWIR band is about 100 ms. If you observe five (5) or six (6) waves in a single SWIR, the AC light source is the cause.

Waves can also occur if the lamp reflector and/or cover glass behave as a white-light interferometer. Solution: remove the glass and/or use a more diffuse reflector.

B.1.11 What are these upward or downward spikes in VNIR data?

These result from artificial light sources, in particular fluorescent lights.

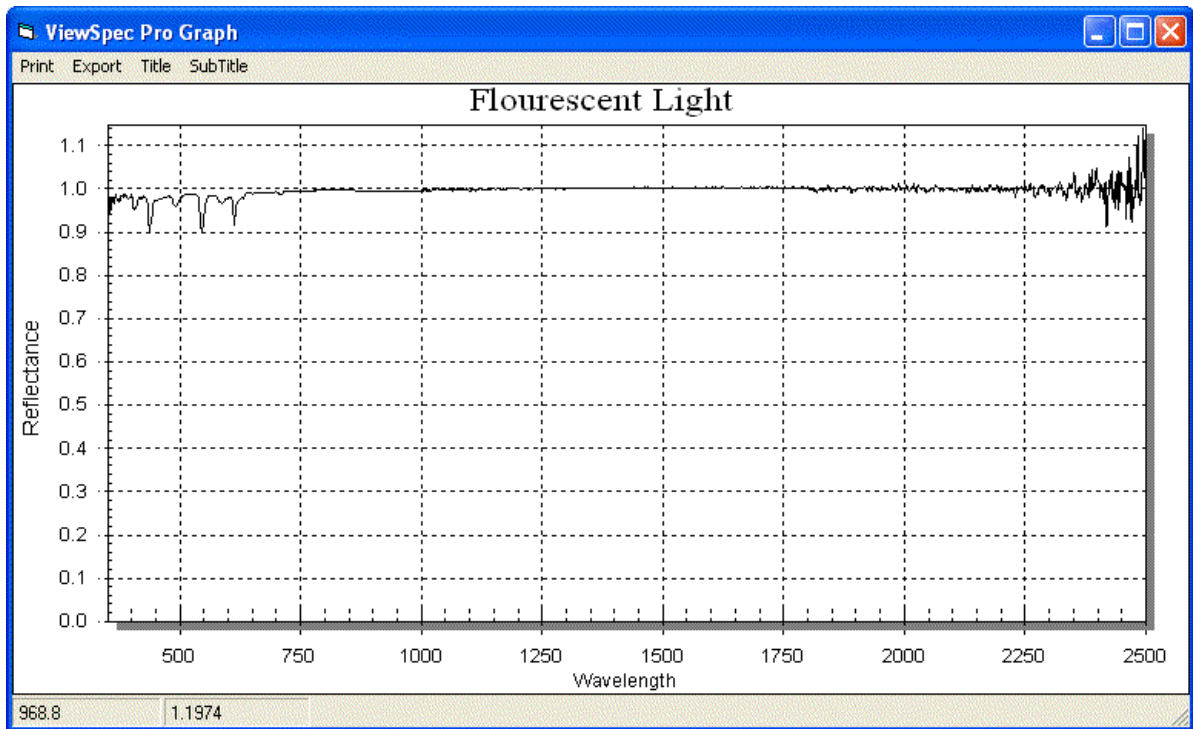


Figure B-1 Spectrum of fluorecent lights

B.1.12 What are these steps in my data?

Light is brought into the instrument by various combinations of the strands in the fiber optic cable. Each strand has its own field-of-view (FOV). When the cable is held close to the sample, each strand views slightly different portions of the sample. This can result in the stepped data.

Stepping of data is common when the foreoptic has a lens, and less common when using the bare optics or sampling devices.

B.1.13 What can cause more noise in my data from last time?

Many factors can cause noise in your data from one session to another. Noise in a measurement is related to the instrument, the signal level, and noise in the light source. Many times the appearance of noise is actually a decrease in the strength of the signal, as opposed to an increase in noise.

Under normal operating conditions, noise visible in a spectrum is always the result of a trade-off between the inherent noise in the system and the signal.

Check how the external fiber optic cable is aligned with the internal cable for changes in signal through-put.

Also, broken fibers in the fiber optic cable can contribute to noise. Perform a fiber optic check to verify.

An increase in noise can be due to a problem in the instrument such as an electronic component malfunction or a grounding problem. This will many times be indicated by a regular pattern to the noise or periodic bursts of noise that are visible over the normal spectra.

B.1.14 Why does the VNIR drop to zero after a Dark Current collection?

When the VNIR drops to zero after a Dark Current has been collected, this may indicate a problem with the VNIR shutter or Dark Current collection routine. To verify, optimize the instrument and listen for the click of the shutter activating. If you do not hear the shutter click, contact technical support and obtain an RMA for sending the instrument in for repairs.

B.1.15 How do I collect a reference with a spot size larger than the Spectralon panel?

Purchase a bigger reference panel or move the sampling accessory closer to it.

If you are indoors using artificial light, make sure that the distance and angle from the reference is the same as will be used for the sample.

B.1.16 How do I set up GPS?

The GPS feature is enabled in the RS³ application software.

To configure the GPS:

Step 1 Select either:

- **GPS->Settings...** pull-down menu item.
- **Alt+G, S** key combination.

RS³ requires the output of the GPS device to be in the National Marine Electronics Association (NMEA) format. Check your GPS device documentation for output and port settings.

Normally, the Baud rate (4800), Data bits (8), Parity (none), and Stop bits (1) settings will remain the same as what is displayed in the window. The **Write to log file** is checked by default. When this field is checked and the GPS fixed data is enabled, an entry will be made to a daily log file (i.e. 050302gps.log) when a spectrum is saved.

- Step 2 In the **GPS Settings** dialog box, enter communication settings appropriate for your instrument controller's port intended for use as well as the GPS device.
- Note: The RS³ settings on the instrument controller and the ports that are used must match the actual port used as well as the settings of the GPS device.
- Step 3 Connect an appropriate cable between the instrument controller's port and the GPS device.
- You may need a serial-USB converter depending upon the instrument controller and GPS device.
- Step 4 Enable the GPS functionality in the RS³ application by selecting either:
- **GPS->Enable...** pull-down menu item.
 - **Alt+G, E** key combination.
- Step 5 Refer to the manual of your actual GPS device for further instructions regarding usage and configuration.

B.1.17 What does a broken fiber mean?

The internal fiber-optic cable is made up of fifty-seven (57) randomly distributed glass fibers.

- 19 fibers are 100 micron for the VNIR region.
- 19 fibers are 200 micron for the SWIR1 region.
- 19 fibers are 200 micron for the SWIR2 region.

Therefore, a single broken fiber in the internal cable results in a ~5% reduction in the signal in a given region.

The external fiber-optic cable is made up of forty (44) 200 micron glass fibers.

- Note: Each broken external fiber results in a few percentage loss of response over the entire range. It is hard to quantify how this affects the response of any given range, because each time the external cable is connected to the spectrometer, its alignment with the fibers of the internal cable will most likely be different.

A few broken fibers is not critical when measuring reflectance, because this is a ratio. Any noise introduced by the broken fibers in measuring the sample were also present when measuring the white reference, and therefore cancels out.

The fibers are protected by a metal spiral inside the black cable casing. If there are kinks in your cable, the fibers are not necessarily damaged. However, if your cable has been crunched so severely that you are able to see

the protective metal spiral, the chances are high that the fibers have been damaged too. Refer to the Fiber Optic Check utility to determine if any fibers are broken.

The fibers can be damaged by coiling the cable up too tightly. If left in a tight coil for longer than a week, the fibers are likely to develop longitudinal fractures that will not be detectable. These fractures in the fiber will cause light leakage, resulting in a weaker signal. The fiber optic cables should be stored by placing them loosely within the netting compartment on the instrument.

WARNING! The fiber cable should never be stored with a bend of less than a 5" diameter for long periods of time.

B.1.18 How long is the battery life?

The charge of the NiMH (Nickel-Metal Hydride) battery lasts about nine (9) hours using a VNIR-only spectrometer with accessory not powered by the instrument in an ambient environment. The charge of the NiMH battery lasts over four (4) hours using a full-range spectrometer with a Contact Probe in an ambient environment.

The Gel Cell rechargeable battery lasts about nine (9) hours using a VNIR-only spectrometer with accessory not powered by the instrument in an ambient environment. The charge of the Gel Cell battery lasts over four (4) hours using a full-range spectrometer with a Contact Probe in an ambient environment.

The Gel Cell rechargeable batteries are rated to give 200 to 300 uses when they are fully drained (discharged to 9 volts) each use. Letting the battery discharge under 9 Volts will shorten the battery's life. If the battery is only used for 50% rated capacity up to 500 uses are possible. In the same way using only 30% capacity can result in up to 1,000 uses.

Note: The limiting factor for how long you can work in the field may be the battery in the instrument controller rather than the spectrometer's batteries.

B.1.19 What type of Ethernet cable can I use for the static IP configuration?

The spectrometer can be plugged directly into the instrument controller using a cross-over, shielded Ethernet cable.

When the spectrometer and instrument controller communicate over a network, use standard shielded Ethernet cables and IP addresses compatible with the network. The spectrometer imposes significant traffic on the network, which can cause packet delays to other users. More importantly, network traffic from other users can negatively impact the reliability of the communication between spectrometer and instrument controller.

B.1.20 Can I install additional software on the instrument controller?

Yes, but with qualifications.

The types of applications that can interfere with the measurement of data are utility programs, network programs, and those working in the background, such as virus checkers.

ASD programs, such as RS³ and Indico, require real-time access to the data that is being streamed from the AgriSpec spectrometer at a high rate of speed. Programs running in the background can cause packets to be lost.

Microsoft Office, image processing programs, and other software applications generally do not interfere with ASD programs, particularly if they are not running and competing for CPU cycles and RAM at the same time that data is being collected from the AgriSpec spectrometer.

B.1.21 Why does the software seem to do unexpected things?

To assure accuracy in the collection and processing of data, the ASD application software is designed to finish its current operation before moving on. The AgriSpec spectrometer outputs a lot of data at a high rate of speed for the ASD application software to collect.

In addition, the ASD application software will *stack up your keystroke entries* and execute them later in the order they were received.

It is important that you wait for the collection to finish before entering the commands to launch another operation. Don't rush into new operations or into issuing new commands until you see the results of the current command.

B.1.22 How long does it take for the AgriSpec spectrometer to warm up?

The warm-up time of the instrument depends on the environment in which it is used.

Only 15 minutes is needed for reflectance measurements, particularly if conserving battery life is desirable.

B.1.23 Where is my serial number?

The serial number is a five (5) digit number located on a label on the rear of the AgriSpec spectrometer where the power switch is. The label also contains the model number and MAC address for the spectrometer.

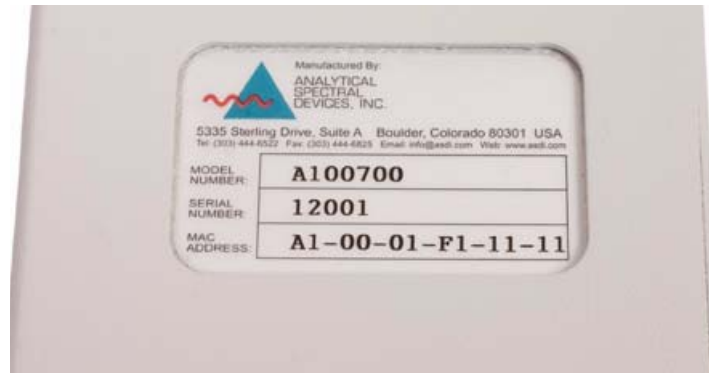


Figure B-2 Label with serial number.

The serial number is also accessible in the RS³ and Indico applications from the splash screen or the **Help->About** menu item.

B.2 Data Conversion

When using the RS³ application, ViewSpec can convert spectral data to ASCII text files.

When using the Indico application, converters are available for Grams/AI, Unscrambler, and ASCII.

Conversion can be done one file at a time. Or, several files can be merged into a single text file, which is a useful feature when inputting data into other analysis programs.

B.3 Gain and Offset

Gains and offsets are determined during optimization by an optimization algorithm in the RS³ application. You can observe this process in the RS³ application by depressing the **OP** button on the top right-hand toolbar. The gains, offsets, and integration time will then appear in the lower left hand corner. These parameters are set by the optimization routine and cannot be set manually.

The **Gain Value** is the inverse of actual gain. A decrease in **Gain Value** will produce a higher gain. The maximum **Gain Value** is 16. Under the low light levels that create such a gain, small changes in **Offset** will translate into large changes in dynamic range.

Offset causes a downward shift (from Y=32,768 DN) in a SWIR spectrum such that the max value of the spectrum and the minimum value of the spectrum (Dark Current - off screen) are equidistant from the line y=32,768. Under ideal conditions, where there's enough light, there will be about 32,768

DN of signal immediately after optimization, +/- 1,000. This means that optimization “creates” about 16,000 DN of dark current and about 16,000 DN of dynamic range. But this dynamic range will disappear with either a slight increase in signal (if the signal started out low to begin with) or a slight increase in dark current (increases with ambient temperature).

B.4 Optimization

The AgriSpec spectrometer must be re-optimized for:

- Every 30 minutes when used with artificial light sources.
- Changes in temperature.

Optimization depends on the response to light in a particular spectral region. A well-optimized instrument will display between 20,000 and 35,000 raw digital numbers, if the source is within the gain and integration time abilities of the system. The magnitude of the VNIR raw DN will be directly dependent on the light intensity.

The Spectralon blank should be used when optimizing and for taking a white reference measurement.

Note: It is important that the position of the reference sample when taking a white reference is as similar as possible to the position for capturing data from the samples.

B.5 Spectrum Averaging (or Sample Count)

If signal levels are low, the only way to increase S/N is by reducing noise through spectrum averaging. However, spectrum averaging takes more time per spectrum.

B.6 Temperature Effects

The AgriSpec spectrometer instruments have thermo-electrically cooled SWIR detectors (to -25 C), significantly reducing noise in those areas.

When the system fails due to extreme temperature, it will fail gradually but the data won't make perfect sense. Shut down and let the instrument cool before operating again.

You may also see the AgriSpec spectrometer saturate in the SWIR 2 region if the instrument is still coming to equilibrium from a cooler environment to a warmer environment. Frequent re-optimizations will address this problem until the instrument comes to equilibrium.

The ASD applications provide a warning when the SWIR detectors are out of thermal regulation.

B.7 Technical Support

If you have any questions or concerns, please contact ASD Inc. by phone, fax, or email:

Phone: 303-444-6522 X-144

Fax: 303-444-6825

email: support@asdi.com

Web: www.asdi.com

ftp: <ftp://ftp.asdi.com>

Technical support is committed to providing you with a timely response to your questions. We will work with you to provide solutions to your applications. Technical support is available to answer your questions Monday thru Friday, 8 am to 5 pm Mountain Standard Time. We will happily respond to your e-mail queries as well.

B.8 W.E.E.E. Compliance



Analytical Spectral Devices supports the W.E.E.E. Regulation. ASD marks the instruments with the symbol at the left to show that this product has entered the market place after August 13, 2005.

If you decide to end the use of the spectrometer, please do not discard it as general waste. ASD will accept the return of your instrument for recycling purposes. Contact ASD Customer Service to arrange for the return of your instrument.

Index

A

access point
 can't connect 46
accessories 5
accessory
 light sources and probes 10
accessory power port 16
AgriSpec 50
array spectrometer 49
ASCII conversion 58

B

baseline collection 51
battery 31
 care 31
battery life 56
broken fiber 55

C

check fiber optic cables 26
cleaning 20
communications 8
computer requirement 6
conversion of data 58

D

dark current in VNIR 54
Dark Current measurement 9
data conversion 58
data sinusoidal wave 52

E

Ethernet 15
Ethernet cable 56
Ethernet interface 5

F

fiber check 26
fiber optic cables 21
fiber optics 21
fiberoptic interface 21
field-of-view 53
FieldSpec 50
Frequently Asked Questions (FAQ) 49

G

gain 58
GPS 54

H

hardware specifications 3

I

illumination 36
instrument controller 6
interface requirement 5

L

LabSpec 50
language settings, Windows 7
LED fiber optic check 26
light sources 19

M

maintenance
 annual 20
materials
 man-made 37

N

near infrared 8
Netcfg 47
NIR 8
noise in data 53

O

offset 58
optimization 51, 59

P

physical specification 3
post-process data 52
power port
 accessory 16
power requirements 3
power supply 13

Q

QualitySpec 50

R

radiance units [52](#)
remote sensing [50](#)
return merchandise authorization [20](#)
RMA [20](#)
rocks [37](#)

S

sample count [51](#)
scanning spectrometer [49](#)
serial number [57](#)
short-wave infrared [8](#)
software requirement [6](#)
software, additional [57](#)
soils [37](#)
spectrograph [49](#)
spectrometer [49](#)
 does not connect [43](#)
 theory of operation [7](#)
spectroradiometer [49](#)
spectrum
 averaging [59](#)
spectrum average [51](#)
steps in data [53](#)
support [ii](#)
SWIR [8](#)

T

technical support [ii](#)
temperature effects [59](#)
TerraSpec [50](#)
trademarks [ii](#)
trigger [17](#)

U

units of radiance [52](#)

V

vegetation [36](#)
ventilation [19](#)
visible/near-infrared [7](#)
VNIR [7](#)
VNIR dark current [54](#)
VNIR spikes [52](#)

W

warm up [57](#)
wavelength specification [4](#)
white reference [9, 39, 51](#)
Windows language settings [7](#)
wireless connection
 can't get [45](#)

loses [47](#)
wireless interface
 disabled [47](#)